


CaseViewer with SlideConverter

Download **Case viewer with slide converter** from the 3DHistech website

Each slide has an MRXS file and a same-name folder (if you rename, change both to the same name)
Right-click on the MRXS file and open with case viewer.

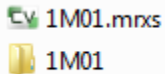

Zoom in and out with control+mousewheel


See the label by using the barcode icon




Select an area to export



The list of annotations should appear automatically, else:

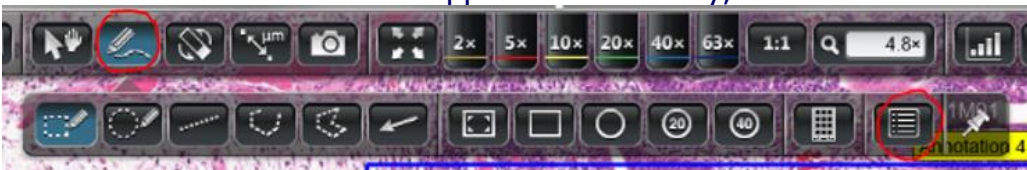


Select an annotation and double click on the "disk" to export

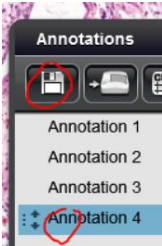


The export dialog should appear.


The icon of the slide converter will appear in the taskbar and may be flashing if the dialog is hidden under a window

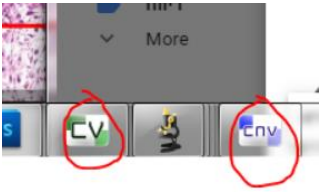

Choose Tiled TIFF

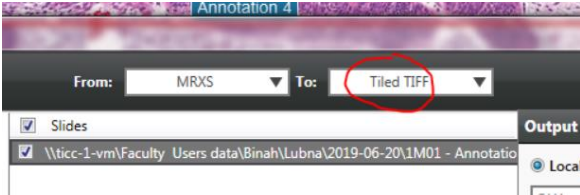

Add slides to queue

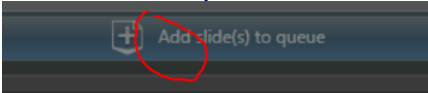

The TIFF file will open with
Windows photo viewer
Photoshop
ImageJ → file → Import → Bioformats (do crop)

In ImageJ you can add scale bar:
Analyze → Tools → Scale bar (do overlay)